



# AQ300

**DISCRETE ANALYZER FOR ENVIRONMENTAL TESTING**

AQ300 is a low cost flexible analyzer that uses the principle of discrete analysis where each test occurs in a separate or discrete reaction vessel.



# AQ300

*AQ300 is ideal for laboratories with smaller workloads requiring varied tests on different samples and when individual results are needed immediately.*

**USEPA, ASTM, ISO**  
and other international regulatory compliant methods are available.

## METHODS INCLUDE

- Alkalinity
- Ammonia
- Chloride
- Cyanides
- Nitrate/Nitrite
- Nitrite
- Phenol
- Phosphate, ortho
- Phosphorus, total
- Silicate
- Sulfate
- Total Kjeldahl Nitrogen

**PLUS MANY MORE**



## HOW DOES THE AQ300 WORK?

The AQ300 robotic sampling arm works in conjunction with a stepper motor-driven syringe that is responsible for aspirating, dispensing and mixing accurate and precise quantities of sample and reagent in miniaturized test tubes called reaction wells.

The sample and reagents are incubated in heated reaction wells until reaction is complete. A single aliquot is then transferred into a 10mm pathlength optical glass cuvette. The absorbance is read on the stationary reactant to ensure the best possible signal to noise ratio.

Each sample is read at the same position in front of the detector in the 10mm glass cuvette, similar to SEAL colorimetric flow systems, known for their high reproducibility and lowest detection limits. This eliminates issues of reaction well variability and scratching found in direct read discrete systems so improves reproducibility. In the SEAL AQ series, liquid is moved not the tray - fewer moving parts equals higher reliability.

Once the absorbance is read, the glass cuvette is thoroughly cleaned and checked, ensuring no carryover or cross contamination.

▶ **As a market leader**, SEAL has over 1,000 applications available and under continual development. Markets include water, wastewater, soil, plant, fertilizer, food and beverage. Please contact us for your specific application.



# AQ300 – The latest in discrete analyzer technology

Designed by chemists for chemists.

## ADVANTAGES AND BENEFITS




- ~ 100% optical glass stop-flow cuvette
- ~ 10 mm optimum path length




- ~ Low cost, disposable wells used for each discrete reaction
- ~ Constant heating and programmable reaction time ensure reaction reaches completion



- ~ Optional integrated cadmium coil reduction for nitrate/nitrite determination
- ~ Cadmium coil is sealed and valve controlled
- ~ In situ coil regeneration is fully automated

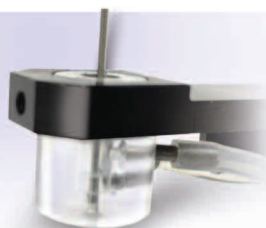

- ~ Unique probe washer for cleaning sample probe between sample and reagent
- ~ Eliminates cross contamination

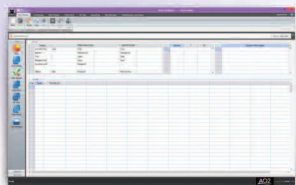

- ~ Highly flexible software designed with user input
- ~ QCPro™ Data Quality System – allows the user to specify QC types, limits and corrective actions



- ▶ **True unattended operation – including ability to run overnight**
- ▶ **Automated standard preparation and dilution of over range samples**
- ▶ **Tests programmable per sample to reduce analysis time**
- ▶ **Add samples after a run has started**
- ▶ **Total volume per test only 500 – 600  $\mu$ L**
- ▶ **LIMS compatible – export in .csv format**
- ▶ **Simple user maintenance**

**SEAL**  
Analytical



PLEASE VISIT

[www.seal-analytical.com](http://www.seal-analytical.com)

or contact your local SEAL Analytical representative about your specific application.

## Colorimetric Nutrient Analyzers

### DISCRETE ANALYZERS



**AQ270**



**AQ300**



**AQ400**

### SEGMENTED FLOW ANALYZERS



**AA100**



**AA500**



**AA3**



**QuAAtro39**

## 50 Years' Experience in Environmental Analysis Built into Every Analyzer

50 years' experience in designing, developing and manufacturing automated wet chemistry analyzers specifically for very low detection levels in environmental applications has helped SEAL to apply the most useful, easy to use features into the SEAL range of Discrete and Segmented Flow analyzers. The SEAL analyzers are widely acknowledged as the best for environmental analysis, giving you everything you need to achieve equal or superior results to the manual and approved laboratory methods the SEAL analyzer replaces.

SEAL Analyzers are monitoring environmental samples in every corner of the globe. They are manufactured in the USA, Germany and the Netherlands. Engineering and chemistry support is provided from SEAL global facilities in USA, Germany, England, the Netherlands and China along with a worldwide network of specialist distributors.

### COMPREHENSIVE SUPPORT

We offer comprehensive applications, technical service and software support.

#### INCLUDING

- ▶ A choice of preventative maintenance and service contracts to meet your specific requirements
- ▶ In-house and online training
- ▶ Guaranteed availability of genuine consumables and spare parts
- ▶ Adaptation of methods to specific requirements such as matrix, range or detection limit
- ▶ Continuous in-house development of software to incorporate new customer requested features

## Robotic Handling Systems

SEAL Robotic Minilab systems for automating sample (pre) treatment in the laboratory – improving your sample handling efficiency. Typical applications include BOD, pH, COD, Alkalinity, and conductivity measurements with options such as decapping/capping, sample splitting, and filtration. Call us about your laboratory needs and we will design a robot to suit you.



**SEAL Minilab 1200**

## Digestion Systems

### FOR METALS AND TKN, TP DIGESTION



**BD50**



**SmartBlock II**



**DEENA 3**

**SEAL Analytical**



[www.seal-analytical.com](http://www.seal-analytical.com)

SEAL Analytical is a global company with offices worldwide - contact us at:

**SEAL Analytical, Inc.**  
6501 West Donges Bay Road  
Mequon, WI 53092  
United States  
Tel: +1 (262) 241 7900  
Fax: +1 (262) 241 7970  
sales@seal-us.com

**SEAL Analytical Ltd.**  
3 Talisman Business Centre  
Duncan Road  
Park Gate  
Southampton, SO31 7GA  
United Kingdom  
Tel: +44 (0) 1489 864 400  
sales.uk@seal-analytical.com

**SEAL Analytical Netherlands**  
ROHASYS BV  
Provincienbaan 4  
5121 DL Rijen  
The Netherlands  
Tel: +31 161 240152  
Fax: +31 161 240153  
info@rohasys.com

**SEAL Analytical GmbH**  
Werkstrasse 5  
D-22844 Norderstedt  
Germany  
Tel: +49 (0)40 60 9292 9-00  
Fax: +49 (0)40 60 9292 9-02  
info.germany@seal-analytical.com

**SEAL Analytical Shanghai**  
Room 413, 12th Building,  
No. 128 Xiangyin Road,  
Shanghai, 200433  
China  
Tel: +86 21 3362 5002  
Fax: +86 21 3362 5002